DI Water in Service Bay

201-DA

To Turn ON DI Water System:

A) OPEN CITY WATER VALVE (YELLOW)

B) OPEN DI WATER VALVES (RED)

C) TURN ON RECIRCULATION PUMP

D) WAIT FOR GREEN RESISTIVITY INDICATOR BEFORE USING DI WATER



City Water Valve (OPEN Position)



DI Water Valves (OPEN Position)

To Turn OFF DI Water System:

A) TURN OFF RECIRCULATION PUMP

B) CLOSE DI WATER VALVES (RED)

C) CLOSE CITY WATER VALVE (YELLOW)

NOTE: THE DI WATER SYSTEM MUST BE SHUT OFF AT THE END OF EVERY WORK/LAB SESSION



Recirculation Pump Switch (ON position)




Resistivity Monitor (Ready Indicated)

DO NOT USE

DI-WATER

IF BOTH WATER TANK

LEDs ARE ORANGE

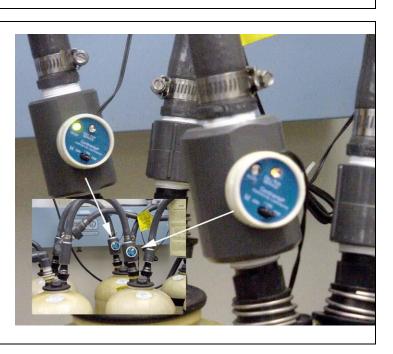

User Manual provided by TJ LaFave, Photos by Cedric Beaulande updated May 9, 2003 (lafatim@iit.edu)